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Foaming suppression system, Platform scale automatic verification system,
Free step sequence

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- Introducing the "foaming suppression system" for handling foamable medical and pharmaceutical products.
- This system is best suited for transporting highly-foamable pharmaceuticals, such as biotechnology-based pharmaceuticals, viscous eye drop solutions, etc.

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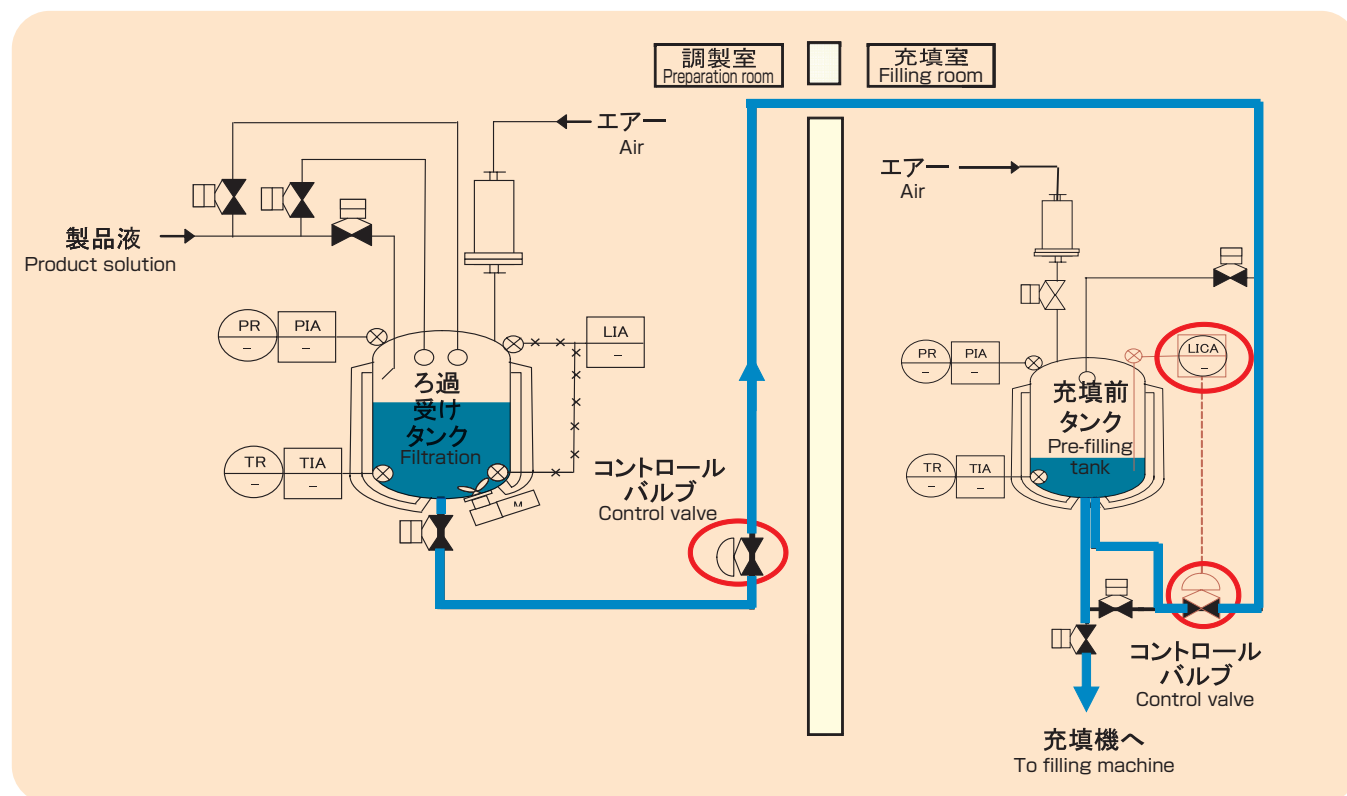
Advantage

- ・Foaming of products during transportation can be suppressed by controlling the degree of opening of two control valves
- ・The pre-filling tank is designed such that a liquid is fed and discharged from the bottom so as to suppress foaming
- ・The filtration receiver tank and the pre-filling tank can be placed on the same floor or on an upper floor and a lower floor.
- ・Highly-accurate pressure control achieves high reproducibility.

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

Platform scale automatic verification system

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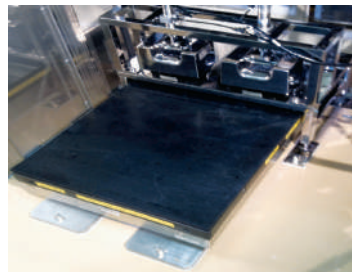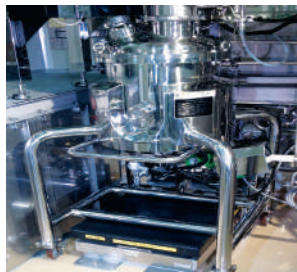Advantage

Avoidance of risks caused by insufficient weighing accuracy

By checking the accuracy of the platform scale before production operation, lot rejection due to insufficient accuracy of the scale can be avoided.

Labor-saving

An actual load for the verification weight is automatically loaded onto the platform scale simply by operating a switch. Heavy manual labor of loading the weight is eliminated. Weighing is carried out by automatically lifting the mobile tank body together with the platform scale.

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Mechanism

- ① A mobile tank is placed on the platform scale.
 - ② The platform scale is lifted by the air cylinder and the mobile tank body is weighed.
 - ③ A verification load weight is automatically loaded by the air cylinder, and accuracy is checked in comparison to the actual load.
- ※ An automatic verification system for stationary tanks is also available.

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- Introducing the "free step sequence" modularization sequence software ideally suitable for multi-product production lines.
- Validation load caused as the result of the addition of production items or change of recipe can be minimized.

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Advantage

- ・ The operation program can be configured by combining the modularized operation program with parameters unique to each item.
- ・ The operation program enables the arbitrary setting of implementation order, combination, operation reservation, etc.
- ・ Operation parameters can be set individually for each item.
- ・ Control change is not necessary when adding a new item. The low development risk CSV category can be applied.

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IWAI PHARMA TECH CO.,LTD.