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- Let us propose the preparation equipment of bonding materials, film coating solutions, etc.

Complete automatic processes of solution preparation from weighing and dissolving to cleaning and sterilizing enable highly-reproducible quality control. This system is optimal for the preparation of a large amount of solution.

移動式調製設備
Mobile preparation equipment

大容量調製設備
Large-capacity preparation equipment

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Advantage

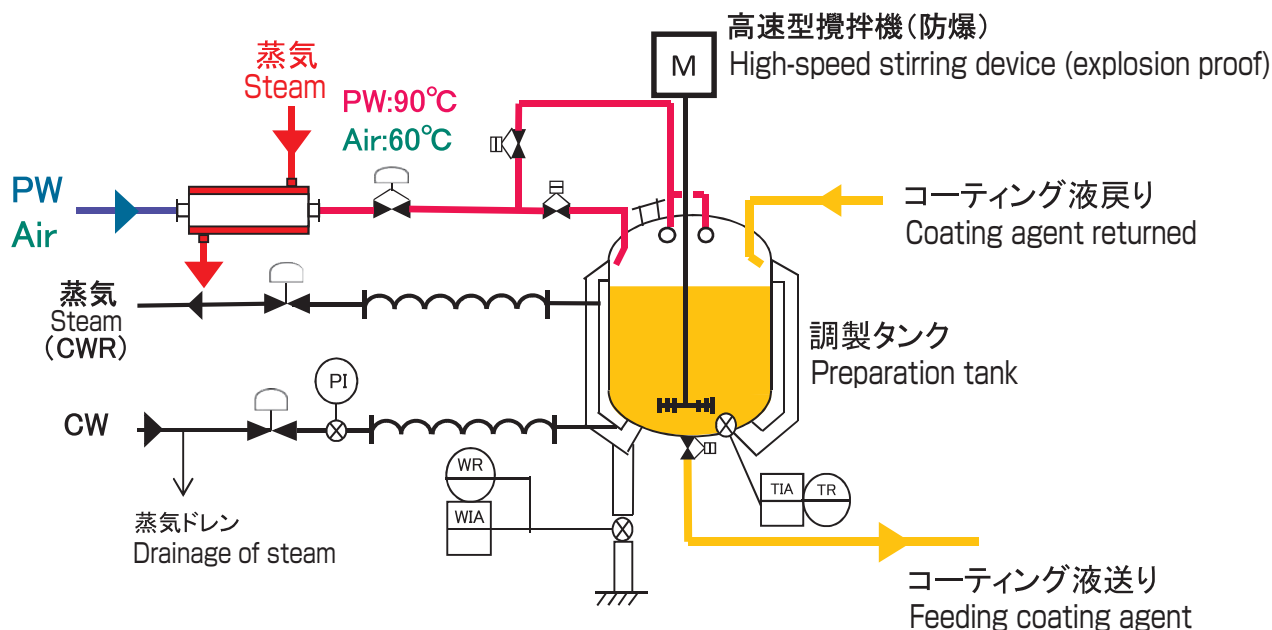
- 1) High-precision weighing
Combination of a load cell and a control valve enables automatically obtaining purified water within the accuracy of ± 0.1 kg.
- 2) Applicable to alcoholic materials and poorly-soluble materials
An explosion-proof stirring device is adopted to ensure dissolution of materials for alcoholic pharmaceuticals.
A high-speed stirring device and special wings are selected for dissolution of poorly-soluble materials to attain dispersion and dissolution.
- 3) Reproducible cleaning, sterilization and drying
Automatic CIP can be conducted with tanks and piping. Warm water sterilization with high-temperature purified water at 95℃. Equipment is dried by thermal air pressure after sterilization to control bio-burden.
- 4) Large capacity
Capable of preparing a large amount of solution up to 2000 L per batch by conducting automatic material weighing, dissolving, heating, cooling, cleaning and sterilizing.

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Solution preparation equipment for solid pharmaceuticals

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Film coating agent preparation equipment flow

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Specifications

Capacity per batch	50 L to 2000 L
Treated items	Binding agent, film coating agent
Type of equipment	Both explosion-proof and non-explosion-proof devices can be used
Accuracy	Accuracy of obtaining water < ± 0.1 Kg Accuracy of temperature regulation < ± 2.0℃
Material of wetted part	Equivalent to SUS304, SUS316L Electrolytically polished (applicable to passivation treatment)
Cleaning, sterilization	Applicable to automatic CIP, applicable to hot water sterilization, applicable to drying

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IWAI PHARMA TECH CO.,LTD.